



Pricelist "BRT inhibitor test for milk producers"

No.	Article	Price
3053	Farm set for beginners	175,00 €
3050	Milk producer test set 25	33,80 €
3055	Milk producer test set 100	121,70 €
3056	Velcro ankle straps, 5 pieces	2,00 €
3057	Digital thermometer	14,29 €
9100	Collie	1,75 €
9112	Parcel delivery	*

* The cost of transportation as well as of insurance will be charged corresponding to expense, except the transport is organised by the consignee.

1 year shelf life of the BRT test tubes, store refrigerated!

AiM – Analytik in Milch Produktions- und Vertriebs-GmbH

Kaiser-Ludwig-Platz 2
80336 Munich
Germany
Phone +49 89-530 75 120
Fax +49 89-530 75 122
info@aimbavaria.com
www.aimbavaria.com



BRT – the inhibitor test

for inspection associations, dairies and milk producers

● fast ●● unambiguous ●●● reliable



Inhibitor free milk?

For sure!



All prices exclusive of VAT and/or taxes. Only the general terms of business of the AIM GmbH are valid. Former pricelists are not effective any more. Subject to modifications and amendments.

This pricelist is valid from January 1, 2012

Analytik in Milch Produktions- und Vertriebs-GmbH
Registration-Court: München HRB 117746. Place of jurisdiction: Munich, Registered office of the company: Munich. Managing directors: Susanne Nüssel and Dr. Christian Baumgartner. Board of directors chairman: Jürgen Ströbel.

Products for Milk Producers





AiM farm set for beginners (Article: 3053)

As a result of the new EU food hygiene regulations milk producers, as food companies, are also subject to quality assurance obligations. They are fully responsible for their products in terms of product liability and preventive consumer protection. With the **farm set for beginners** from AiM GmbH, milk producers can carry out the officially prescribed inhibitor test (Brilliant Black Reduction Test, BRT) themselves on the farm.

Everything you need for the test is contained in a handy bag. You can now test milk from treated cows after the withdrawal time has elapsed as well as milk to be delivered before it is collected by the tanker fast and reliably to ensure that it is free of inhibitors.



Composition:

- Thermo-block for 5 BRT test tubes
- Milk producer test set
- Product information
- Timer
- Edding
- Sample protocol
- Bottle for milk samples
- Bag
- Velcro ankle straps

AiM milk producer test set (Article: 3050/3055)

The milk producer test set may be reordered at any time. It is available in two packing units. Article 3050 contains 25 BRT test tubes, 25 disposable pipettes and the Inhibitor Free Milk (negative control). Article 3055 contains 100 BRT test tubes, 100 disposable pipettes and the Inhibitor Free Milk.



Accessories


Velcro ankle straps to mark treated cows (article 3056), digital thermometer to check the temperature of your thermo-block (article 3057)



Performing the test is easy:

1. Preparation

- Switch the thermo-block on
- Mix sample well
- Mark the test tubes

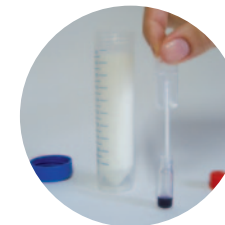

2. Preparing the negative control

Dip the pipette into the milk and activate it by pressing the upper air chamber. Pipette the negative control (Inhibitor Free Milk) into a test tube and then close the tube with the cap.



3. Preparing the samples

Important: Use a different pipette and BRT test tube for each sample.

4. Incubate the test tubes

65 °C

Incubate the test tubes until the test medium in the tube with the negative control changes from blue to yellow.

2:30 hours +/- 15 minutes

5. Interpretation of the results

- left = Negative control
(Inhibitor free sample)
- middle = Suspicious sample
- right = Sample contains inhibitors

Milk with positive/suspicious results must not be delivered!

