



## Pricelist "BRT inhibitor test for milk producers"

No.	Article	Price
3053	Farm set for beginners 25	182,00 €
3054	Farm set for beginners 48	196,00 €
3050	Milk producer test set 25	35,00 €
3058	Milk producer test set 48	49,00 €
3055	Milk producer test set 100	126,00 €
3056	Velcro ankle straps, 2 pieces	2,50 €
3057	Digital thermometer	17,90 €
3059	Timer	19,90 €
9100	Collie	1,80 €
9112	Parcel delivery	*

### AiM – Analytik in Milch Produktions- und Vertriebs-GmbH

Kaiser-Ludwig-Platz 2  
80336 Munich  
Germany  
Phone +49 89-530 75 120  
Fax +49 89-530 75 122  
info@aimbavaria.com  
[www.aimbavaria.com](http://www.aimbavaria.com)



## BRT – the inhibitor test

for inspection associations, dairies and milk producers

● fast ●● unambiguous ●●● reliable



Inhibitor free milk?

**For sure!**



\* The cost of transportation as well as of insurance will be charged corresponding to expense, except the transport is organised by the consignee.

**1 year shelf life of the BRT test tubes, store refrigerated!**


All prices exclusive of VAT and/or taxes. Only the general terms of business of the AiM GmbH are valid. Former pricelists are not effective any more. Subject to modifications and amendments.

This pricelist is valid from January 1, 2014

Analytik in Milch Produktions- und Vertriebs-GmbH  
Registration-Court: München HRB 117746. Place of jurisdiction: Munich, Registered office of the company: Munich. Managing directors: Susanne Nüssel and Dr. Christian Baumgartner. Board of directors chairman: Günther Felßner.

Products for Milk Producers



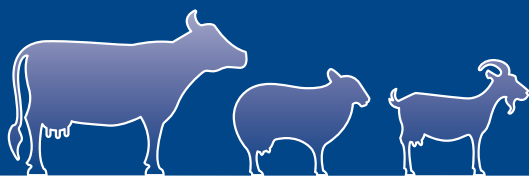

## AiM farm set for beginners (Article: 3053/3054)

As a result of the new EU food hygiene regulations milk producers, as food companies, are also subject to quality assurance obligations. They are fully responsible for their products in terms of product liability and preventive consumer protection. With the **farm set for beginners** from AiM GmbH, milk producers can carry out the officially prescribed inhibitor test (Brilliant Black Reduction Test, BRT) themselves on the farm.

Everything you need for the test is contained in a handy bag. You can now test milk from treated cows after the withdrawal time has elapsed as well as milk to be delivered before it is collected by the tanker fast and reliably to ensure that it is free of inhibitors.



### Composition:

- Thermo-block for 5 BRT test tubes
- Milk producer test set 25 resp. 48
- Product information
- Timer
- Edding
- Sample protocol
- Bottle for milk samples
- Bag
- Velcro ankle straps

## AiM milk producer test set (Article: 3050/3055/3058)

The milk producer test set may be reordered at any time. It is available in three packing units. Article 3050 and 3055 contain 25 resp. 100 glass BRT test tubes with screw caps, article 3058 contains 48 plastic BRT test tubes. All sets contain a sample protocol, the corresponding number of disposable pipettes and the Inhibitor Free Milk (negative control).



### Accessories


Velcro ankle straps to mark treated cows (article 3056), digital thermometer to check the temperature of your thermo-block (article 3057), timer to check incubation time of the BRT test tubes (article 3059)



## Performing the test is easy:

### 1. Preparation

- Switch the thermo-block on
- Mix sample well
- Mark the test tubes

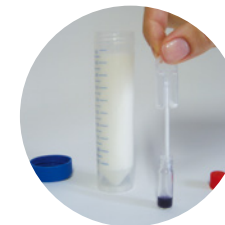

### 2. Preparing the negative control

Dip the pipette into the milk and activate it by pressing the upper air chamber. Pipette the negative control (Inhibitor Free Milk) into a test tube and then close the tube with the cap.

### 3. Preparing the samples

Important: Use a different pipette and BRT test tube for each sample.



### 4. Incubate the test tubes

65 °C

Incubate the test tubes until the test medium in the tube with the negative control changes from blue to yellow.

2:30 hours +/- 15 minutes

### 5. Interpretation of the results

left = Negative control  
(Inhibitor free sample)

middle = Suspicious sample

right = Sample contains inhibitors

Milk with positive/suspicious results must not be delivered!

