



BRT test tubes – this is the way to get an unambiguous and reliable result

Inhibitor testing starts with sampling.
The BRT Inhibitor Test produces correct results only with high-quality milk samples which were taken properly.



Sampling

- Mix whole milking well before sampling
- Avoid contamination
- Use clean sample containers
- Take sample with a clean ladle



Quality of samples:

- Only test fresh samples
- Ensure that all samples are kept refrigerated
- Avoid heat and direct sunlight
- Don't test sour samples



Performing the test

- Store the test tubes in the refrigerator
- Mark the samples and the test tubes
- Switch on the heating block about 10 mins in advance
- Mix the milk sample and don't take samples from the cream
- Use disposable pipettes and bin them after use to avoid contamination
- Use a negative control
- Pay attention to the incubation time (about 2 ½ hrs)
- Record the results in the sample protocol to ensure traceability



Farm tests are for your own use and are no official tests or results.